

**IMCO Corporation Ltd AB**

PO Box 21195  
SE -100 31 Stockholm  
Sweden  
[www.imcocorp.se](http://www.imcocorp.se)

## PRODUCT SPECIFICATION

<b>Product No.</b>	<b>FkTRXR-03*FS</b>
<b>Product description:</b>	<b>Fluorescent substrate for Thioredoxin reductase assay with FkTRXR-03*</b>
<b>Lot No.</b>	<b>01</b>
<b>Specifications:</b>	<b>Fluorescent substrate</b> Lyophilized fluorescent substrate <i>Preparation procedure:</i> Dissolve the content in 1.0 ml distilled water. Store at -20°C
<b>Reference:</b>	See methods description for FkTRXR-03* in <a href="http://www.imcocorp.se">www.imcocorp.se</a>
<b>Storage:</b>	Keep at -20°C
<b>Examples of use:</b>	A substrate for measurement of thioredoxin reductase using FkTRXR-03* kit from IMCO. See <a href="http://www.imcocorp.se">www.imcocorp.se</a>

**NOTE!** Make sure to reconstitute all of the lyophilized material in the ampoule. After adding water, close the screw-cap again, shake vigorously, then centrifuge shortly, in order to recover liquid from tube cap and walls.